

Instructions for use

2x qPCR Probe BLUE MasterMix (high ROX)

1 ml, 100 reactions

Optimized 2x PCR MasterMix for all probe-based qPCR assays; including ROX reference dye (high concentration), blue coloured

1. Description

Our **2x qPCR Probe BLUE Master Mix** is an optimized ready-to-use mixture for probe-based assays such as TaqMan®, Beacons and MGBs. It contains a modified fast HotStart Taq DNA Polymerase, dNTPs and MgCl₂ combined in an optimized buffer system for Real-Time PCR / qPCR applications except primers, probe and template DNA / cDNA. The Master Mix can be used not only for expression analysis but also for genotyping, copy-number analyses and all sorts of probe-based quantitative PCRs. The **2x qPCR Probe BLUE MasterMix** contains ROX reference dye at a concentration of 1 µM (in its final concentration 500nM) and a blue dye for convenient pipetting.

For research use only. Not approved for use in clinical or *in vitro* diagnostics.

2. Applications

2x qPCR Probe BLUE MasterMix is recommended for use in all probe-based Real-Time PCR / qPCR applications such as TaqMan®, Beacons and MGBs. The optimized buffer system provides fast kinetics and highly efficient target amplification with low DNA template amounts and even for difficult templates. The **2x qPCR Probe BLUE MasterMix** contains all components, you just need to add primers, probe and template DNA / cDNA. The **2x qPCR Probe BLUE MasterMix** is not only suitable for expression analysis, but also for genotyping, copy-number analysis and all sorts of probe-based quantitative PCR assays. Using our **2x qPCR Probe BLUE MasterMix** not only reduces contamination risks, but is also time saving, highly reproducible and very easy to prepare.

3. Set contents

2x qPCR Probe BLUE MasterMix in **2x reaction buffer** containing the modified HotStart Taq polymerase, 0.4 mM each dNTP and an optimized buffer system containing MgCl₂, 1 µM ROX and a blue dye.

Reagent	Amount	Lid colour
2x qPCR Probe BLUE MasterMix (100 reactions)	1 tube, 1 ml	brown

Note: The **2x qPCR Probe BLUE MasterMix** is also available without or with low concentrations of ROX.

4. Protocol

Before you start: Thaw the tube and invert the MasterMix 5-6 times to ensure mixing of the solution. Do not vortex! After thawing spin the tube briefly!

Note: Reactions can be conveniently set up at room temperature.

Recommended reaction mixture per well:

Components	Apply for 20 μ l reaction	Apply for 10 μ l reaction	Final concentration
2x qPCR Probe BLUE MM	10 μ l	5 μ l	1x
Forward primer	variable (e.g. 2 μ l)	variable (e.g. 1 μ l)	0.1 – 0.4 μ M
Reverse primer	variable (e.g. 2 μ l)	variable (e.g. 1 μ l)	0.1 – 0.4 μ M
Probe	variable (e.g. 2 μ l)	variable (e.g. 1 μ l)	0.2 – 0.4 μ M
Template DNA	variable	variable	1 pg – 10 ng/reaction
Sterile dest. water	adjust to 20 μl	adjust to 10 μl	

- **3-Step PCR protocol**

Step	Time	Temperature
Initial denaturation	2 minutes	92-95 °C
30 - 45 cycles		
Denaturation	5 seconds	92-95 °C
Annealing	5 seconds	60 °C (depends on primers)
Extension	5 – 20 seconds	72 °C

- **2-Step PCR protocol**

Step	Time	Temperature
Initial denaturation	2 minutes	92 - 95 °C
30 – 45 cycles		
Denaturation	5 seconds	92 - 95 °C
Annealing/Extension combined	5- 30 seconds	60 °C (depends on primers)

Note: For maximum efficiency and specificity, annealing temperatures, as well as extension time, primer/probe concentration and template DNA concentration may need to be optimized.

5. Storage conditions

Long-term storage: at -20°C (stable for about 24 months).

Short-term storage: at 4°C (stable for about 3 months).

However, short term storage for a few hours or up to 3 days at room temperature will not affect the performance of the MasterMix.

Product is not covered by pending or issued patents or may have certain limitations. To our best knowledge, this product does not provide any conflict with pending or issued patents.